

# GUINEA PIG SNUGGLE

## INSTRUCTIONS FOR USE



### Part of the SNUGGLE Family of Comfortable Restrainers

Congratulations on the purchase of your GUINEA PIG SNUGGLE.  
Please make sure you have purchased the correct size for your animal.

You are now ready to start.

1. First open your SNUGGLE on a flat surface with the rear extension (1) facing you.
2. Place the animal on the SNUGGLE.
3. Close by pulling tail flaps (1) up over the back, (male (hook) Velcro will be facing upwards).
4. Fold body flap (2) over tail flaps to secure the animal in the SNUGGLE.
5. Fold body flap (3) to close the restrainer and to adjust for individual fit.
6. Finally close flaps (4) at the neck to secure the animal.

The SNUGGLE may be used in many configurations for various testing and sampling procedures.

All SNUGGLES are protected by US patent # 5,975,028. All rights are reserved.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to [info@lomir.com](mailto:info@lomir.com).

### Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC  
Canada J7V 7M4  
Tel: 514-425-3604  
Toll Free: 1-877-425-3604  
Fax: 514-425-3605  
Email: [info@lomir.com](mailto:info@lomir.com)

### Lomir Biomedical Inc.

C/O Agenda Life Sciences  
UK PO Box 24 Hull  
HU128YK, UK  
Tel: 084566060556  
Email: [info@lomirbiomedical.co.uk](mailto:info@lomirbiomedical.co.uk)

### Lomir Inc.

PO Box 778  
213 West Main Street  
Malone, NY  
USA 12953  
Tel: 518-483-7697  
Toll Free: 1-877-425-3604  
Email: [info@lomir.com](mailto:info@lomir.com)



# CARE INSTRUCTIONS

## FOR YOUR SNUGGLES

To clean SNUGGLES, start by soaking for up to an hour in a quaternary ammonium disinfectant, this removes heavy soiling, any odor as well as disinfecting. Do not use chlorine bleach. Then proceed to wash using industrial detergent (such as cage wash detergent) or laundry soap in a washing machine.

Similar results can be achieved by placing SNUGGLES in a cage washer. All disinfectants and detergents are used at the manufacturers' recommended concentrations.

Temperature – Best results will be achieved in warm water.

For best results, close any Velcro fasteners, this protects the Velcro from pulling or attaching to other fabrics and SNUGGLES. SNUGGLES may be machine dried.

All SNUGGLES are protected by US patent # 5,975,028. All rights are reserved.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to [info@lomir.com](mailto:info@lomir.com).

### Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC  
Canada J7V 7M4  
Tel: 514-425-3604  
Toll Free: 1-877-425-3604  
Fax: 514-425-3605  
Email: [info@lomir.com](mailto:info@lomir.com)

### Lomir Biomedical Inc.

C/O Agenda Life Sciences  
UK PO Box 24 Hull  
HU128YK, UK  
Tel: 084566060556  
Email: [info@lomirbiomedical.co.uk](mailto:info@lomirbiomedical.co.uk)

### Lomir Inc.

PO Box 778  
213 West Main Street  
Malone, NY  
USA 12953  
Tel: 518-483-7697  
Toll Free: 1-877-425-3604  
Email: [info@lomir.com](mailto:info@lomir.com)