

# JACKETS CARE

## Instructions



To clean jackets, start by soaking for up to an hour in a quaternary ammonium disinfectant, this removes heavy soiling, any odor as well as disinfecting. Do not use chlorine bleach. Then proceed to wash using industrial detergent (such as cage wash detergent) or laundry soap in a washing machine. Similar results can be achieved by placing jackets in a cage washer. All disinfectants and detergents are used at the manufacturers' recommended concentrations. Temperature – Best results will be achieved in warm water. For best results, close any Velcro fasteners, this protects the Velcro from pulling or attaching to other fabrics and jackets. Make sure that Spandex inserts or jackets incorporating Spandex are not washed in hot water. Jackets may be machine dried.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to [info@lomir.com](mailto:info@lomir.com).

### Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC  
Canada J7V 7M4  
Tel: 514-425-3604  
Toll Free: 1-877-425-3604  
Fax: 514-425-3605  
Email: [info@lomir.com](mailto:info@lomir.com)

### Lomir Biomedical Inc.

C/O Agenda Life Sciences  
UK PO Box 24 Hull  
HU128YK, UK  
Tel: 084566060556  
Email: [info@lomirbiomedical.co.uk](mailto:info@lomirbiomedical.co.uk)

### Lomir Inc.

PO Box 778  
213 West Main Street  
Malone, NY  
USA 12953  
Tel: 518-483-7697  
Toll Free: 1-877-425-3604  
Email: [info@lomir.com](mailto:info@lomir.com)