



BUNNY SNUGGLE

INSTRUCTIONS FOR USE

Part of the SNUGGLE Family of Comfortable Restraints

Congratulations on the purchase of your new BUNNY SNUGGLE the safest form of rabbit restraint. Please assure you have purchased the correct size for your rabbits.

You are now ready to start your first stress free rabbit procedure.

1. First open your SNUGGLE on a flat surface with the rear extension (1) facing you.
2. With your subjects head tucked under your arm place the subject on the SNUGGLE facing (8).
3. While holding the subject firmly, raise panel (1) and hold with the thumb and forefinger of your hand.
4. Close flap (2) over (1) then proceed to cover (2) with (3). Assure a firm closure.
5. Now fasten (5) over (4) behind the neck. Next, cover the eyes with flap (7) secured over flap (6).
This will keep the subject calm during procedure.

Panels (7) and (4) as well as (6) and (5) may be interfaced for a more secure hold of the head (gavage, eye movement etc.)

The SNUGGLE may be used in many configurations for various testing and sampling procedures.

All SNUGGLES are protected by US patent # 5,975,028. All rights are reserved.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7V 7M4
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 084566060556
Email: info@lomirbiomedical.co.uk

Lomir Inc.

PO Box 778
213 West Main Street
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com



CARE INSTRUCTIONS

FOR YOUR SNUGGLES

To clean SNUGGLES, start by soaking for up to an hour in a quaternary ammonium disinfectant, this removes heavy soiling, any odor as well as disinfecting. Do not use chlorine bleach. Then proceed to wash using industrial detergent (such as cage wash detergent) or laundry soap in a washing machine.

Similar results can be achieved by placing SNUGGLES in a cage washer. All disinfectants and detergents are used at the manufacturers' recommended concentrations.

Temperature – Best results will be achieved in warm water.

For best results, close any Velcro fasteners, this protects the Velcro from pulling or attaching to other fabrics and SNUGGLES. SNUGGLES may be machine dried.

All SNUGGLES are protected by US patent # 5,975,028. All rights are reserved.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7V 7M4
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 084566060556
Email: info@lomirbiomedical.co.uk

Lomir Inc.

PO Box 778
213 West Main Street
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com