

DOG SLING

SUIT & FRAME



FRAME

Put the base on the table with the small holes facing up. Take the up-rights and line them up with the holes on the base plate with corresponding holes on the base. From underside, insert the stainless bolts (1/4" -20/ 1 1/2" stainless bolts) through the holes. Use the wing nuts to tighten the bolts. Repeat for all 16 bolts. Once both sets of up-rights are on, you may tighten using a 7/16" open wrench. Pass the cross-member of the frame through the zipper sleeve of the sling suit. Repeat for the other side. Place the cross-member in the up-rights and tighten them by passing 1 1/2" bolts through the corresponding holes using the wing nuts to tighten the bolts.

SLING SUIT

Introduce the animal to the sling. Detach the body of the suit from the zipper sleeve. Undo the 2 right hand zippers to the jacket. Place the animal in the jacket and close the zippers. Zip the animal in the sling suit to the zipper sleeves on the frame.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7V 7M4
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 084566060556
Email: info@lomirbiomedical.co.uk

Lomir Inc.

PO Box 778
213 West Main Street
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com