

LARGE ANIMAL

FRAME ASSEMBLY INSTRUCTIONS

WHEELS

Insert the wheel assembly into the frame leg to the desired height. Secure the assembly in place using the 1 1/2" bolt and wing nut. Hand-tighten. It is recommended that the two locking wheels be located at opposite corners of the frame. (Figure 1)

HANDLE ASSEMBLY

Insert stud screws in threaded holes provided. (Figure 2-3)
Put handles in between screws. (Figure 4-5)

TIE DOWNS

Insert single end or ropes into cleats. (Figure 6)



Figure 1



Figure 2-3



Figure 4-5



Figure 6

The fabric sling or hammock is fitted to the frame by passing the side poles through the channels on either long side of the hammock. It is recommended that the animals be acclimated to the sling in advance of any procedures.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7V 7M4
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 084566060556
Email: info@lomirbiomedical.co.uk

Lomir Inc.

PO Box 778
213 West Main Street
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com