



P.E.T 2027

CARE & USAGE

PRIMATE ENRICHMENT

Thank you for your purchase of the 2027 P.E.T. This enrichment device is designed to be used with New World Primates. The device is manufactured from food grade polypropylene; all fasteners are stainless steel. The principal of operation is that the food is located inside the device and is accessed by the subject turning an elliptical wheel and aligning the holes.

CLEANING:

The P.E.T. environmental enrichment device may be cleaned by a number of techniques. For example, emersion in a bleach or disinfectant solution followed by rinsing, placing in a cage or tunnel washer. They may be sterilized using an autoclave.

FILLING:

For cleaning and filling the 2027 P.E.T., unscrew the knob on top of the unit and remove the large disk. The device is filled with selected food items and the disk and upper knob are both replaced.

FOOD/TREATS:

The 2027 device can be filled with a range of food items. The following are suggestions: dried fruits, fruit sauce or jellies, nut butter, propriety treats or diet (e.g. Lab Diet's Callitrichid diets 5LK6 or 5LK7). Mealworms could also be used with the outer disks preventing the mealworms from escaping.

PLACEMENT OF DEVICE:

The 2027 can be hung inside an enclosure or cage.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7V 7M4
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 084566060556
Email: info@lomirbiomedical.co.uk

Lomir Inc.

PO Box 778
213 West Main Street
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com