



# P.E.T 2029

---

## CARE & USAGE

### PRIMATE ENRICHMENT

Thank you for your purchase of the 2029 P.E.T. This enrichment device is designed to be used with New World Primates. The device is manufactured from food grade polypropylene, all fasteners are stainless steel.

#### CLEANING:

The P.E.T. environmental enrichment device may be cleaned by a number of techniques. For example, emersion in a bleach or disinfectant solution followed by rinsing, placing in a cage or tunnel washer. They may be sterilized using an autoclave.

#### FILLING:

For filling and cleaning the 2029 P.E.T., unscrew the front and remove the disk(s). For sauce, jelly and gels use a large spatula to quickly fill all cavities.

#### FOOD/TREATS

The 2029 device can be filled with a range of food items. The following are suggestions: dried fruits, fruit sauce or jellies, nut butter, propriety treats or diet (e.g. Lab Diet's Callitrichid diets 5LK6 or 5LK7). Clear H<sub>2</sub>O gels such as LabGel® could also be used.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to [info@lomir.com](mailto:info@lomir.com).

### Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC  
Canada J7V 7M4  
Tel: 514-425-3604  
Toll Free: 1-877-425-3604  
Fax: 514-425-3605  
Email: [info@lomir.com](mailto:info@lomir.com)

### Lomir Biomedical Inc.

C/O Agenda Life Sciences  
UK PO Box 24 Hull  
HU128YK, UK  
Tel: 084566060556  
Email: [info@lomirbiomedical.co.uk](mailto:info@lomirbiomedical.co.uk)

### Lomir Inc.

PO Box 778  
213 West Main Street  
Malone, NY  
USA 12953  
Tel: 518-483-7697  
Toll Free: 1-877-425-3604  
Email: [info@lomir.com](mailto:info@lomir.com)



# P.E.T 2029

## CARE & USAGE

### PLACEMENT OF DEVICE:

The 2029 can be hung inside an enclosure or cage or can be mounted on the front of a cage. All devices are supplied with two stainless steel eye bolts for mounting, an additional set of two threaded holes at the base of the device allow for additional attachments.

The 2029 device is supplied with two 2" diameter plastic disks. You will note that these disks are two different thicknesses; the correct disk is selected according to how you are using the device and whether you are using one or both rotating plates:

- ½" x 2" disk is to be use with a single plate.
- ¼" x 2" disk is to be used when both plates are in use.

To dismantle for cleaning, changing the plates and filling with food/treats, hold stainless hex bolt with a wrench and unscrew plastic disk on the outer portion of the device.

Animals may use various tools to forage in the smaller holes located in the top and bottom trays and when the plate with the smaller holes is used.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to [info@lomir.com](mailto:info@lomir.com).

### Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC  
Canada J7V 7M4  
Tel: 514-425-3604  
Toll Free: 1-877-425-3604  
Fax: 514-425-3605  
Email: [info@lomir.com](mailto:info@lomir.com)

### Lomir Biomedical Inc.

C/O Agenda Life Sciences  
UK PO Box 24 Hull  
HU128YK, UK  
Tel: 084566060556  
Email: [info@lomirbiomedical.co.uk](mailto:info@lomirbiomedical.co.uk)

### Lomir Inc.

PO Box 778  
213 West Main Street  
Malone, NY  
USA 12953  
Tel: 518-483-7697  
Toll Free: 1-877-425-3604  
Email: [info@lomir.com](mailto:info@lomir.com)