



# P.E.T 2031

## CARE & USAGE

### PRIMATE ENRICHMENT

Thank you for your purchase of the 2031 P.E.T. This enrichment device is designed to be used with New World Primates. The device is manufactured from food grade polypropylene, all fasteners are stainless steel.

#### CLEANING:

The P.E.T. environmental enrichment device may be cleaned by a number of techniques. For example, emersion in a bleach or disinfectant solution followed by rinsing, placing in a cage or tunnel washer. They may be sterilized using an autoclave.

#### FILLING:

For filling and cleaning the 2031 P.E.T., unscrew the front and remove the disk. Use a tongue depressor to fill cavities. For sauce, jelly and gels use a large spatula to quickly fill all cavities.

#### FOOD/TREATS:

The 2031 device can be filled with a range of food items. The following are suggestions: dried fruits, fruit sauce or jellies, nut butter, propriety treats or diet (e.g. Lab Diet's Callitrichid diets 5LK6 or 5LK7). Clear H2O gels such as LabGel® could also be used.

#### PLACEMENT OF DEVICE:

The 2031 can be hung inside an enclosure or cage or can be mounted on the front of a cage. All devices are supplied with two stainless steel eye bolts for mounting, an additional set of two threaded holes at the base of the device allow for additional attachments.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to [info@lomir.com](mailto:info@lomir.com).

### Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC  
Canada J7V 7M4  
Tel: 514-425-3604  
Toll Free: 1-877-425-3604  
Fax: 514-425-3605  
Email: [info@lomir.com](mailto:info@lomir.com)

### Lomir Biomedical Inc.

C/O Agenda Life Sciences  
UK PO Box 24 Hull  
HU128YK, UK  
Tel: 084566060556  
Email: [info@lomirbiomedical.co.uk](mailto:info@lomirbiomedical.co.uk)

### Lomir Inc.

PO Box 778  
213 West Main Street  
Malone, NY  
USA 12953  
Tel: 518-483-7697  
Toll Free: 1-877-425-3604  
Email: [info@lomir.com](mailto:info@lomir.com)