

RODENT INFUSION

Rodent Infusion Set

- Rodent jacket
- Rodent tether
- Rodent swivel
- Swivel tether connector
- Swivel retainer



IMPORTANT

A test run using saline on all assembled equipment should be performed prior to your experiment to assure proper flow and a system free of leaks or blockages

ASSEMBLY DIRECTIONS

Before commencing, ensure that the jacket is the correct size for the animal. The tether length is selected according to the dimensions of the cage in use. Affix swivel holder (5) to cage using appropriate model. Diagram shows model RS 01 that clips to the outside of the cage. Insert swivel (4) into swivel holder (5), tighten thumb-screw (6), **DO NOT OVER TIGHTEN**. Attach connector block (3) to the swivel, (4) tightening the Allen screw with the Allen key provided. At this stage, pass the catheter from the animal through the tether (2) and fix the tether to the jacket (1) using the Velcrotab. Hooks provided on the jacket are there to aid in jacket adjustment. They should be tied with dental floss or surgical wire. For additional security, it is extremely important that this connection in the system does not turn. Affix the tether (2) into the connecting block (3), tightening the Allen screw with the Allen key provided. Attach the catheter to the swivel (4), using an over under fashion to create a loop in the system.

Lomir Biomedical Inc.

95 Rue, Notre-Dame-de-l'Île-Perrot, QC
Canada J7V 7M4
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.


C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 08456600556
Email: info@lomirbiomedical.co.uk

Lomir Inc.

PO Box 778
213 West Main Street
Malone, NY
USA 12953
Tel: 518-483-7897
Toll Free: 1-877-425-3604
Email: info@lomir.com

RODENT INFUSION

Rodent Infusion Set

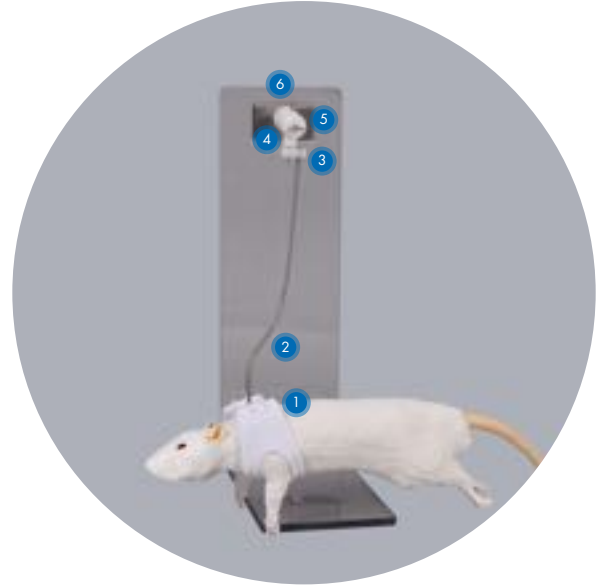

ROUTINE MAINTENANCE

On a regular basis, check all connections in the system. This may be undertaken at the same time as routine animal observations are made.

STERILIZATION

Preferred method of sterilization is ethylene oxide, or “wet” immersion sterilization. It is not recommended to autoclave swivels.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7V 7M4
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 084566060556
Email: info@lomirbiomedical.co.uk

Lomir Inc.

PO Box 778
213 West Main Street
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com