



STERILIZATION OF LOMIR SWIVELS AND OTHER EQUIPMENT

Recommended as first choice for sterilization is Ethylene oxide (ETO) and similar proprietary methods of gas sterilization. You may also use “wet” -immersion sterilization using a range of disinfectant solutions. All solutions would be prepared and used as per the manufacturers’ recommendations. It is **NOT** recommended to autoclave or add high temperatures to any of the above procedures. Lomir swivels should never be sterilized by irradiation.



Should you have any questions or need more information please contact our Customer Service Department toll free in North America at 1-877-425-3604 or send an e-mail to info@lomir.com.

Lomir Biomedical Inc.

95 Huot, Notre-Dame-de-l'Île-Perrot, QC
Canada J7V 7M4
Tel: 514-425-3604
Toll Free: 1-877-425-3604
Fax: 514-425-3605
Email: info@lomir.com

Lomir Biomedical Inc.

C/O Agenda Life Sciences
UK PO Box 24 Hull
HU128YK, UK
Tel: 084566060556
Email: info@lomirbiomedical.co.uk

Lomir Inc.

PO Box 778
213 West Main Street
Malone, NY
USA 12953
Tel: 518-483-7697
Toll Free: 1-877-425-3604
Email: info@lomir.com